

EyeSpray First-Aid Eye and Face Wash

Model 850.1003

1] INSTALLATION INSTRUCTIONS

- Determine a visible location, close to the danger.
- Install the rack on the wall (screws not supplied), so that the spray bottles are at a height of 1.1 m, as for any eyewash, which corresponds to a distance of 1.15 m between the bottom of the closed rack and the floor.
- Install the bottles according to replacement instructions below, making sure the nozzles are oriented upwards.
- Close the cover of the rack, leaving the red flag inside. Fix the red sealing across the cover and the baseplate so that it will necessarily be broken when activating the eyewash. Also fix the label indicating the expiry date (3 years after installation).

Regularly inform and train the users and the security team as to the protocols to follow in case of emergency.



2] REPLACEMENT OF THE SPRAY BOTTLES

After each use, new bottles must be installed, and the seal must be replaced. The red flag indicates that the eyewash has been activated and that new bottles must be installed.

- Open the cover of the rack, however not completely, leaving a space - using a pen or an object - thus allowing the spray bottles not to be activated.
- Dismount the old bottles from their casing and remove them from the rack.
- Make sure to have two new bottles available, as well as a white label to indicate the expiry date and a red sealing sticker.
- Write down the expiry date on the white label. The bottles are to be replaced three years after installation.
- Remove the cap from the new bottles. Install the new bottles in the rack, making sure the nozzles are facing upwards and the bottles are firmly fixed in their casing. A supplied sticker reminds about the need to correctly orient the nozzles.
- Close the cover of the rack, leaving the red flag inside, thus showing the rack has not been opened after replacement of the bottles.
- Fix the red sealing across the cover and the baseplate so that it will necessarily be broken when activating the eyewash. Also fix the white label with the expiry date, preferably on the side of the rack.

A periodical control is necessary to make sure that the seal is not broken, that the red flag is not visible, and that the expiry date has not been reached.

These instructions can be found, and a video can be seen on www.eyespray.ch



SCAN FOR
INSTRUCTION VIDEOS

LaboSafe SA	www.eyespray.ch - www.labosafe.ch	GIA-Premix www.giapremix.ch
Rue de Saint-Nicolas 10 CH-2000 Neuchâtel	EyeSpray : Installation and Maintenance Instructions	
Phone: +41 32 756 96 96	info@labosafe.ch	En / 2022-11